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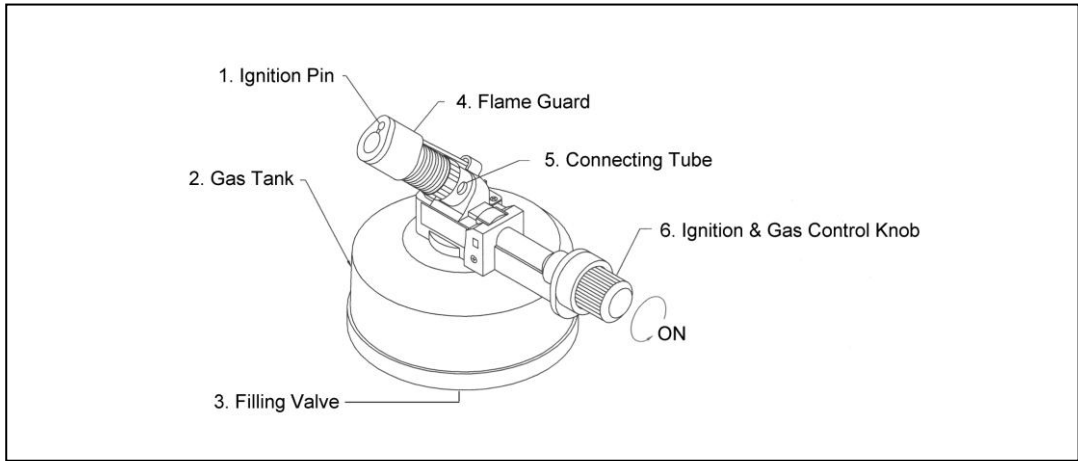
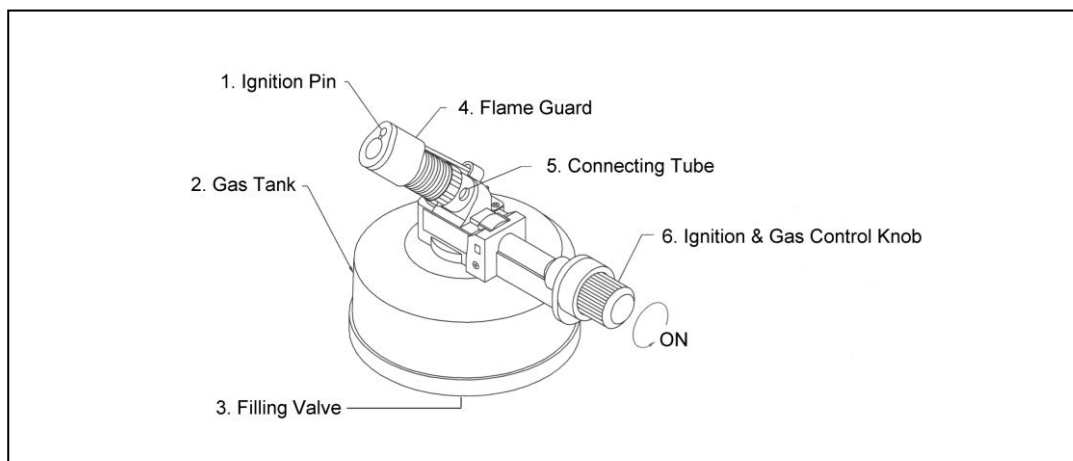
Dragon 220 TILT HEAD LAB BURNER

ATTENTION: This unit does not contain any butane gas when purchased. Please follow below to make it operational.

1. FUEL FILLING AND REFILLING

- a) Make sure the gas burner is switched off before filling.
- b) Reverse the unit upside down,
- c) Insert the container's nozzle vertically in a downward motion into the FILLING VALVE ③ on the bottom of the unit.
- d) When the fuel begins to overflow, remove the nozzle from the filling valve to stop the filling operation. It means the unit is full of gas.
- e) Allow a few minutes for the fuel to stabilize in the fuel tank.

2. IGNITION

- a) Please adjust the position of Air Control Sleeve to align the small hole on the sleeve with the hole on the CONNECTING TUBE. ⑤
- b) Turn the IGNITION BUTTON ⑥ to the left (counterclockwise) about 90 degree.
- c) Press the IGNITION BUTTON ⑥ to light the flame.

3. SHUT-OFF

Turn the IGNITION BUTTON ⑥ to right (clockwise) to decrease until the end to shut off the flame.

4. FLAME LENGTH ADJUSTMENT

After the ignition, adjust the CONNECTING TUBE ⑤ to adjust the flame shape.

5. WARNING

- a) Do not touch FLAME GUARD ④ when hot.
- b) Keep out of reach of children.
- c) Do not drop. Do not operate or store in temperature exceeding 50°C.
- d) Do not try to disassemble the unit.